MEDICLAVE

4047725

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GENERAL INFORMATION

- 1) *Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributor immediately.*
- 2) *Do not install or use the equipment without reading this handbook before.*
- 3) *This handbook must always be attached to the equipment and it must be available for all users.*
- 4) *If you have any doubts or enquiries, please contact with your supplier or J.P. Selecta's technical service.*
- 5) **IMPORTANT! J.P. SELECTA WILL NOT ACCEPT ANY APPARATUS TO BE REPAIRED IF IT IS NOT DULY CLEANED.**
- 6) *If any modification, elimination or lacking in maintenance of any device of the equipment by the user transgress the directive 89/655/CEE, the manufacturer is not responsible for the damage that can occur.*
- 7) *Do not use the apparatus with liquids which can give off vapours capable of making explosive mixtures.*

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CONTENTS

GENERAL INFORMATION
CONTENTS
PACKING LISTS
ACCESSORIES
TECHNICAL FEATURES
EQUIPMENT DESCRIPTION
INSTALLATION
OPERATION
MAINTENANCE
SPARE PARTS
ELECTRICAL DIAGRAM
GUARANTEE
"EC" CONFORMITY DECLARATION

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The standard equipment consist of the following components:

Descripción / Description	Código / Code	Cant. / Qty.
MEDICLAVE	4047725	1
Cestillo / Basket	1000596	1
Manual de instrucciones / Instruction Manual	80xxx	1

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* Baskets in perforated plate of AISI 316 stainless steel:

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Voltage supply 115-230V 50/60 Hz according to the characteristics plate indications.

Código / Code	4047725	
Capacidad litros / Capacity litres	12	
Medidas útiles cm. Chamber dimensiones cm.	Ø	24.5
	Alto / Depth	25
Medidas exteriores cm. Overall dimensions cm	Alto / Height	50
	Ancho / Length	38
	Fondo / Width	50
Consumo / Consumption W	1000	
Peso / Weight Kg	18	

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EQUIPMENT DESCRIPTION

For sterilisation temperatures of 121°C and 134°C (of and 2 Kg/cm²) fulfilling the current standards for pressure apparatus, certified and signed by the industry services. Top and outer casing in AISI 304 stainless steel, chamber, lid and lock system in AISI 316 stainless steel.

Great safety operation, with automatic air exhaust at start for sterilization cycle. Built in a safety valve and a safety thermostat for excess temperature or lack of water.

INSTALLATION

Place the autoclave near a suitable electric mains.

Level the autoclave with the legs.

Always use decalcified water.

Fix a hose in the water/steam exit nipple of the machine, and also fix the other end to a drain or deposit without **OBSTRUCTING THE FLOW**.

Do not install the autoclave in places where inflammable liquids are stored or in zones with special protection.

Before opening the cover **ALWAYS** check that there is **NOT PRESSURE**. The pressure gauge indicator installed in the control panel must read 0kg/cm².

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Be sure that the voltage supply is the same as the one indicated on the characteristics plate.

Do not use the autoclave if it is not earthed.

NEVER exceed the maximum pressure of 2 Kg/cm²

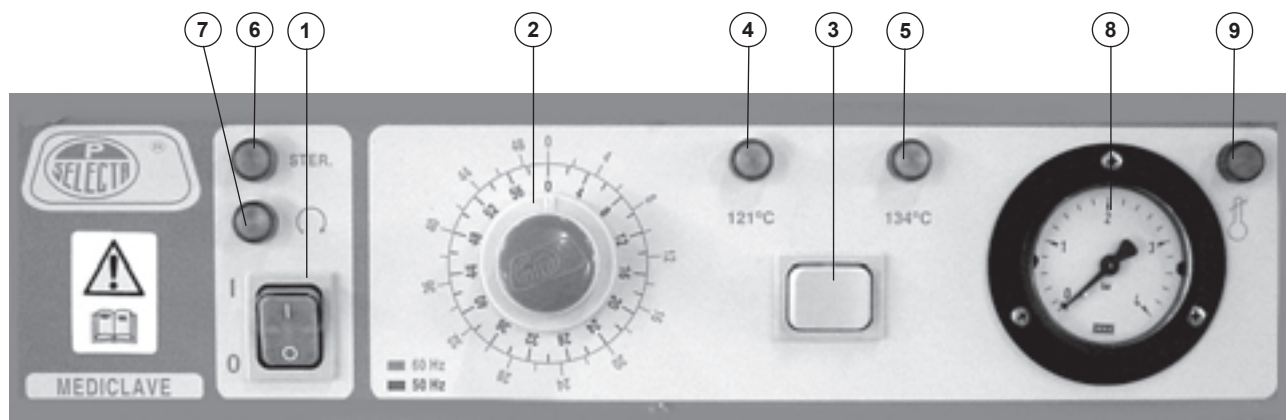
From time to time revise the functioning of the safety valve.

If you change the plug bear in mind the following:

Blue cable: Neutral.
Brown cable: Phase.
Yellow/green cable: Earth.

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STARTING UP:

1. Check that the STEAM valve (10) and the DRAIN-AGE (11) valve are closed.
2. Fill the tank of the autoclave (better with decalcified water) up to the tube rack level.
3. Place the material to be sterilized.
4. Close the cover (12) pressing strongly.
5. Select the sterilization temperature with the selector knob (3). At the moment of sterilisation the relation between temperature and pressure of the saturated steam will be the following: (See table 1).
6. Select the sterilization time pressing the switch knob (2).
7. Press the main switch POWER (1), at the same time the yellow indicator lamp (4) or (5) will light up depending of the selected temperature. The autoclave has an automatic air exit valve which remains open during heating up to 110°C, allowing air removing from the autoclave, and on reaching the previously mentioned temperature, it automatically closes.
8. Once the set temperature is reached the light (6) becomes ON, and the preset time starts count. At the end of preset time light (7) comes ON and autoclave stops operation.

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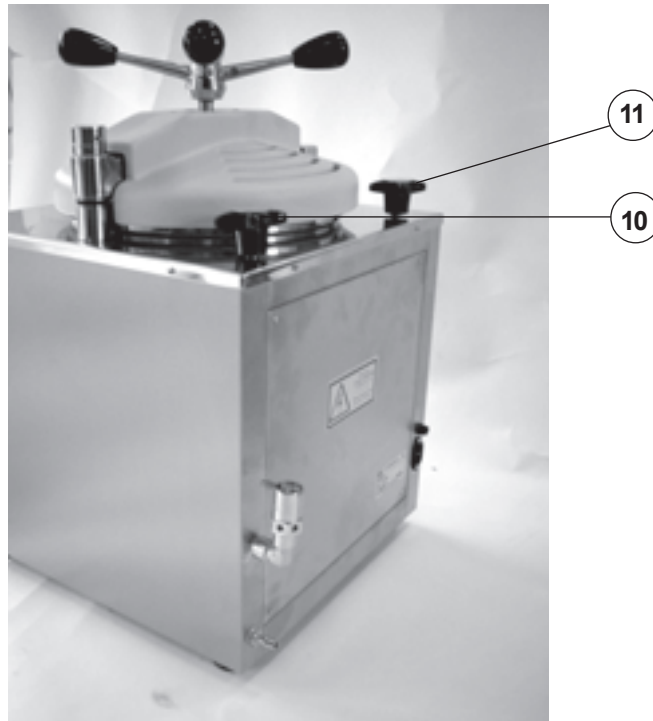
9. Stops autoclave with (1).

ATTENTION!!! DO NOT OPEN THE COVER OF
THE AUTOCLAVE UNTIL THE PRESSURE
GAUGE IS AT «0» AND THE STEAM VALVE IS
OPEN

10. To empty the inner water, place the handle (11) to
DRAINAGE position.

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11. If SOLIDS (gauzes, gloves, cottonwool, instruments, etc...) have been sterilized, steam can be cleared quickly by opening the STEAM valve (10) and emptying the water from the tank with the DRAINAGE valve (11).

12. If LIQUIDS (liquid medicines, closed culture medium, bottles, etc...) have been sterilized, the pressure must return to the ambient temperature alone or in extreme case by VERY CAREFULLY opening the STEAM valve (10).

13. The RAPID DECOMPRESSION OF LIQUIDS provokes the breakage of flasks, spilling of liquids, etc..., with the resulting obstruction of tubes in the lower part of the autoclave.

14. In any case, SOLIDS or LIQUIDS, before open the lid, check that manometer (8) indicates 0 bar and open the valve (10) VAPOR.

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When abnormal increase of the temperature or a considerable decrease in the water level occurs (with the consequent overheating of the heater element), the safety thermostat starts working blocking the heating of the machine and the ALARM indicator lamp (9) lights up. To deblock the machine proceed as follows:

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1. *Disconnect the autoclave from the mains.*
2. *Open the STEAM valve.*
3. *Let it cool.*
4. *Open and fill the tank with water up to the working level.*
5. *Reset the safety thermostat (12) placed at the lower front part.*
6. *Start up the autoclave.*

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If the problem continues, you must call the Technical Service.

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Said valve is gauged to initiate the opening at 2.2kg/cm², and only begins to work if the pressure switch fails or by an error selecting it.

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MAINTENANCE

Before removing the cover disconnect the apparatus from the mains.

For safety reasons do NOT modify the wiring system. The manipulation of the internal control elements of the apparatus by by unauthorized personnel can cause irreparable damage.

CLEANING:

For the cleaning of the different parts use the following products:

Cleaning of stainless steel: Alcohol

Cleaning of plastic: Alcohol with cotton duster.

Cleaning of pipes with lime: Solution of Hydrochloric acid 50% and water 50%.

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Annual revision of the whole function of the autoclave by qualified personnel is advisable.

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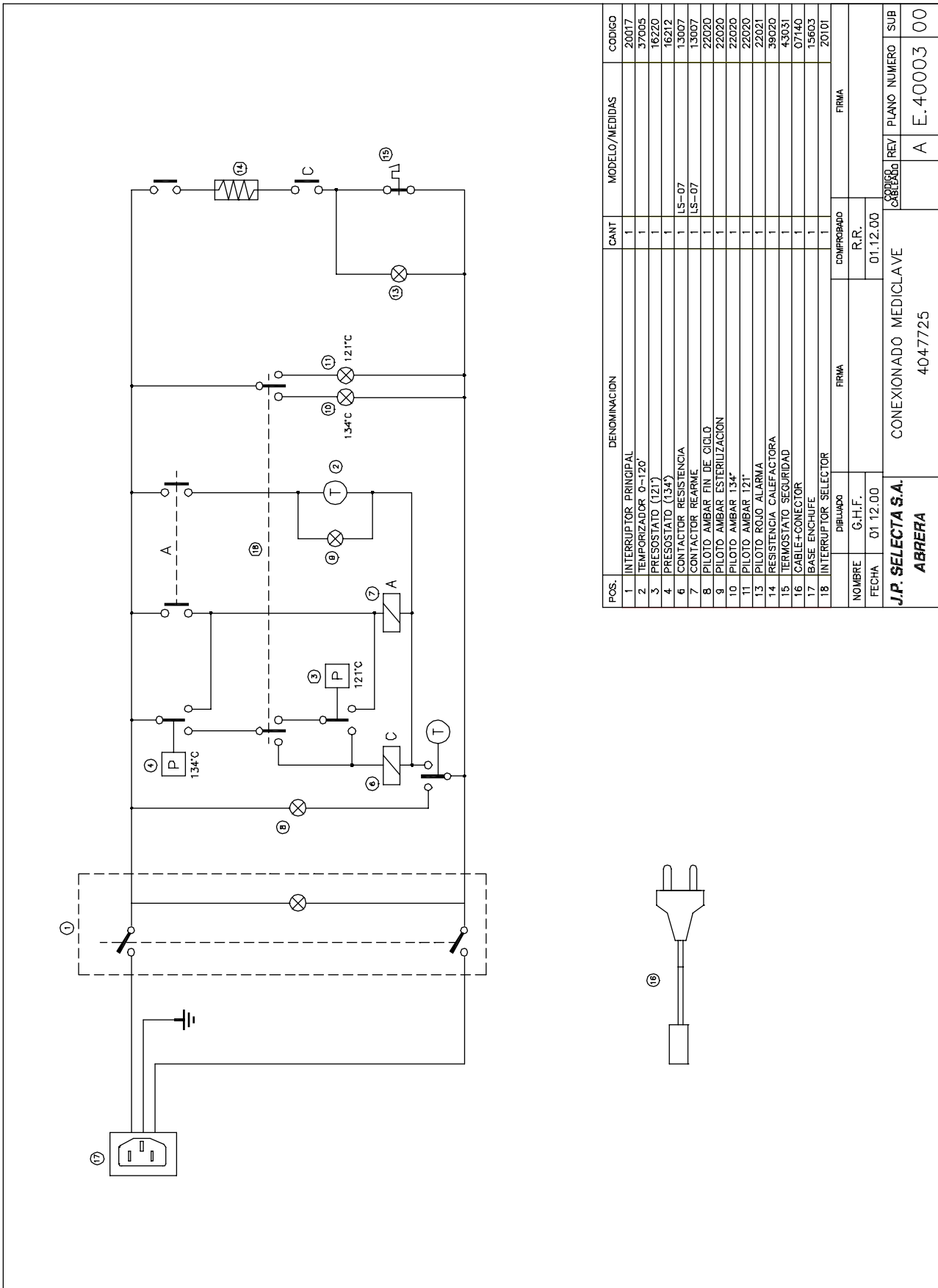
SPARE PARTS

To guarantee the safety of the equipment, the spare parts must be purchased from J.P.SELECTA,S.A.

Descripción / Description	4047725
Gradilla / Tube rack	5602
Manómetro / Manometer	16002
Presostato / Pressure switch	16212
Presostato / Pressure switch	16220
Interruptor principal / Main switch	20017
Interruptor selector / Selector switch	20061
Junta tapa / Gasket cover	21059
Piloto verde / Greene neon lamp	22017
Piloto ámbar / Ambar neon lamp	22020
Piloto rojo / Red neon lamp	22021
Válvula seguridad / Safety valve	34200
Reloj / Timer	37013
Resistencia / Heating element	39159

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GUARANTEE

This product is guaranteed for one year. The guarantee does not cover damage caused by incorrect use or causes beyond the control of J.P. SELECTA, S.A.

Any manipulation of the apparatus by personnel not authorized by J.P. SELECTA, S.A. automatically cancels the guarantee.

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