



## Paraffin dispenser 43900



### Instruction manual

1. Fill paraffin container with paraffin and close lid - 12 - .
2. Apparatus supply with voltage.
3. Turn temperature pre-selecting knob - 11 - to position 10 (90°C).
4. Turn on master switch - 3 -. Both pilot lamps - 4 - + - 5 - are lighting.  
Due to the thermostatic temperature control the yellow pilot lamp - 5 - will temporarily go out and light up again.
5. After melting of the paraffin (about 1 hour) turn down to a lower temperature depending upon the sort of paraffin used (for example, 62°C - turn temperature pre-selecting knob - 11 - between position 8 + 9).  
The selected temperature will be maintained by the thermostatic temperature control.
6. The hose from your vacuum pump must be connected to the hole - 13 -, in order for the gases to be syphomed out.
7. In order to discharge the desired quantity of paraffin, depress operating lever of the filling tap - 14 - Quantity to be discharged can be regulated by finger pressure.
8. Turn off the apparatus only by using the master switch - 3 - .

**This equipment should not used in explosion-endangered areas.**

MEDAX GmbH & Co. KG  
Kieler Str. 211, D-24768 Rendsburg  
Tel.: +49 4331-136-870, Fax: +49 4331-136-838  
e-mail: [info@MEDAX-Kiel.de](mailto:info@MEDAX-Kiel.de)  
[www.MEDAX-Kiel.de](http://www.MEDAX-Kiel.de)

## Spare Part List

Pos.	Description	Part-No.	
1	Power plug with integrated fuse	MEN 28100	
2	Fuse, 2,5 A	MEN 11700	F1 / F2
3	Tumbler switch	MEN 11000	S1
4	Pilot lamp, green	MEN 10600	H1
5	Pilot lamp, yellow	MEN 01100	H2
6	Fuse of temperature	MEN 01700	F3
7	Condenser	MEN 25400	C1
8	Thermostat	MEN 00100	B1 / B2
9	Bottom heating element, 300W	MEN 10000	E1
10	Tap-heating, 50W	MEN 05600	E2
11	Knob	MEN 01800	

## Electrical Diagram

