




Paraffin Tissue Floating Bath, round model 24900, 25900

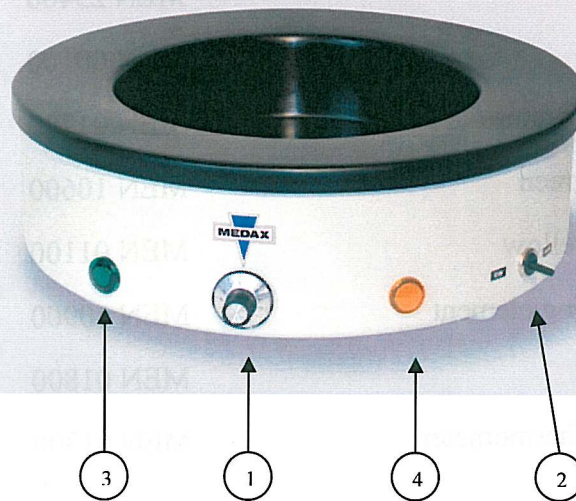

Instruction manual

1. Before using, fill up the container with distilled water.
2. Fix the holder with thermometer on the edge of the water bath.
3. Connect the power cable.
4. Adjust the control knob "1", lower or higher temperatures can be adjusted by turning the control knob counterclockwise or clockwise.
5. Turn on master switch "2". Both pilot lamps "3" and "4" are lighting. The temporary flashing of lamp "4" is subject to the regulator.
The temperature of the water can be displayed on the thermometer.
6. **Make sure, that you have always enough distilled water in the bath.**
7. The Floating Bath should only switch off by using the power-switch "2".

Attention


Never heat the apparatus with empty bath!

This equipment should not used in explosion-endangered areas.

Spare Part List

Pos.	Description	Part-No.	
1	Power plug with integrated fuse	MEN 28100	
2	Fuse, 2,5 A	MEN 11700	F1 / F2
3	Swith	MEN 11000	S1
4	Condenser	MEN 25400	C1
5	Thermostat	MEN 00100	B1
6	Fuse of temperature	MEN 01700	F3
7	Pilot lamp, green	MEN 10600	H1
8	Pilot lamp, yellow	MEN 01100	H2
9	Bottom heating element	MEN 10900	E1
10	Knob	MEN 01800	
11	Mechanical thermometer	MEN 01300	

Electrical Diagram

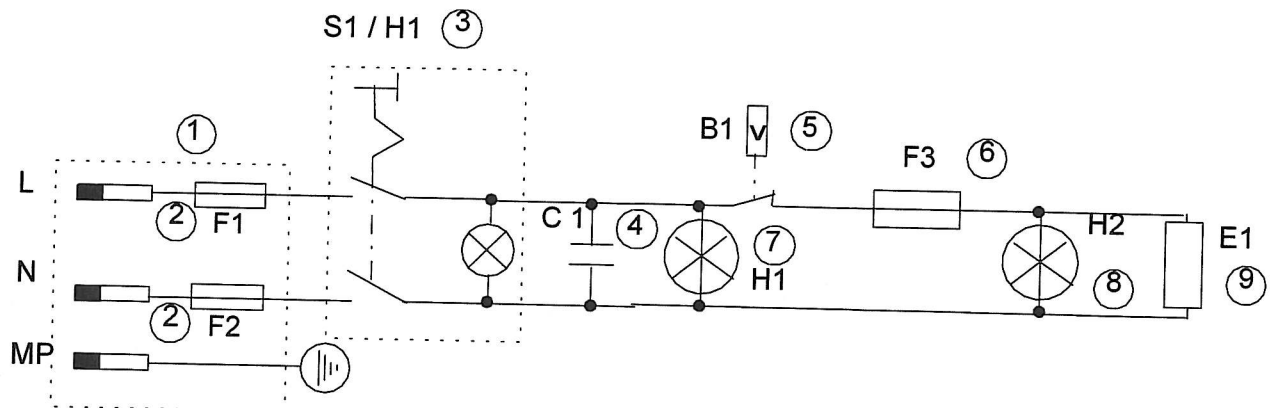

MEDAX GmbH & Co. KG

Kieler Str. 211, D-24768 Rendsburg, Germany

Tel.: +49-(0)4331-136-870, Fax: +49-(0)4331-136-838

e-mail: info@MEDAX-Kiel.de

www.MEDAX-Kiel.de

Subject to technical change, Dec. 05